

How to interface JAGA16 with your electrodes

Jan, 2017

How to interface JAGA16 with your electrodes

- JAGA16 device ends with femal Millmax 18 pins (MILL-MAX 853-93-100-10-001000)
- Below we show several ways to attach your electrodes
- See also JAGA16 Millmax-pin-orientations.pdf on our website.



Millmax EIB


- You can add 4 x tetrodes to this Millmax EIB



Millmax EIB

Millmax-Omnetics adatper


- If your electrodes end already with Omnetics

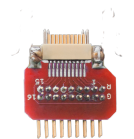

Milmax Omnetics adapter

Millmax Male Pins

- Quick method. You can solder your wires into this Millmax male pins

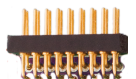

Milmax male pins